



Diode Laser for Nanosecond Pulses with Programmable Shapes

- Center wavelength: 1062 nm
- Pulse width between 1 ns and 330 ns with programmable pulse shape
- Repetition rates up to 1 MHz
- External trigger input
- Various operation modes: bursts, sequences, burst sequences
- Synchronization output
- Signaling for fiber amplifiers



Applications

- Seeding of fiber amplifiers
- Material processing

preliminary preliminary preliminary

Diode Laser for Nanosecond Pulses with Programmable Shapes

The PPL 400 is a stand-alone, computer controlled laser system for the generation of nanosecond pulses with programmable pulse shapes. It is connected to the host computer via a standard USB connection.

The central part of the laser system are up to 4 individual programmable waveform generators. Each generator outputs a combination of a flat top square pulse and a sawtooth with adjustable amplitude and slew rate. Start and stop times of the pulse can be chosen within one of three time bases in each channel. The outputs from all waveform generators are combined, and the sum signal is used to directly modulate the output of a laser diode. The optical output power of the laser is fully voltage controlled which allows a much better extinction ratio than electrooptically modulated sources.

The PPL 400 can be programmed through an easy to use Windows[™] control software. The software provides a graphical preview of the output of all waveform generators as well as the combined signal that drives the laser diode and thus corresponds to the optical pulse's temporal modulation. Last settings are saved inside the PPL 400 to allow stand-alone operation.

The PPL 400 consists of a mainframe with a power supply, an oscillator module (SOM 828), one or two dual channel waveform generation modules (SWM 828) and a voltage controlled laser module (VCL 828) with polarization maintaining fiber output.



Waveform generation module - SWM 828

Channel 1
pulse amplitude [‰]
210 0 500 1000
pulse start [%]
290 0 500 1000
ramp slew (%)
0 500 1000
ramp start [‰]
290 0 500 1000
wave stop [%)
900 0 500 1000
timebase 330 ns

The PPL 400 can be equipped with one or two waveform generation modules (SWM 828). Each module includes two channels with six variable parameters for the generated waveform:

- the time base (33 ns, 100 ns or 330 ns)
- the start time and the stop time of the pulse within the chosen time base
- the amplitude of the flat top square pulse
- the start time and ramp slew of the sawtooth within the chosen time base

The combination of these parameters allows generation of a large variety of pulse shapes, from the very basic flat top square pulse to highly sophisticated multipeak patterns.

The output of each waveform generator is connected to one input of the voltage controlled laser module (VCL 828).

The SWM 828 also features a "status control" sync output which can be connected to, e.g., an amplifier system, in order to indicate proper operation. The SWM 828 can work at any repetition rate up to 1 MHz (depending on selected time base).

Voltage controlled laser module - VCL 828

The VCL 828 module is used to generate the laser emission based on the driving signals from the SWM 828 module. The output from one or two SWM 828 modules is internally combined and used to directly control the output power of the laser emission, which is highly linear with respect to the input voltage within a certain range. The laser has a lasing threshold of typically 1 mW and a maximum output power of typically 100 mW. Within this range the laser power corresponds to the control voltage between 10 mV and 1 V. At 0 V input voltage the laser is completely switched off and thus provides a very high extinction ratio between the pulses.

preliminary preliminary preliminary

Oscillator module - SOM 828

The oscillator module of the PPL 400 is identical to the oscillator module of the PDL 828 "Sepia II" from PicoQuant.

It has eight (individual) trigger outputs (channel) that can be addressed individually. For example, channels can be combined to be activated at the same time or each channel can be individually activated in a sequence. To enable burst operation, multiple pulses can be output from one channel (or from combined channels) before the next channel becomes active.

Working principle

Generally speaking, the oscillator module works in a rotary fashion, which means that the programmed sequence of channel 1 through channel 8 must be completed before channel 1 becomes active again as illustrated in figure 1.

All channels can therefore be activated individually in a sequence. Adjacent channels can also be combined for simultaneous operation. The channels can, however, not be combined in any arbitrary fashion as the rotary working principle must be preserved.

For the PPL 400, only two channels of the SOM 828 are used for providing trigger signals to the waveform generators. The six other channels are thus free and can for example be used to define pauses between two bursts

Repetition rate

The oscillator module provides a wide range of internal, user selectable repetition rates between 196 kHz and 80 MHz, which are derived from an internal crystal oscillator (80, 64 or 50 MHz) along with a frequency divider (1 to 255). The SOM 828 can also be triggered using external signals at any repetition rate up to 80 MHz. The frequency divider is also active on external trigger signals, which therefore permits to realize virtually any repetition rate between single shot and 80 MHz.

The highest truly usable repetition rate for the SOM 828 within a PPL 400 device is, however, defined by the chosen time base on the waveform generation module and therefore limited to 30 MHz. At higher repetition rates, the pulses are no longer separated and the laser will switch to continuous wave operation.

Defining bursts and combining channels


The oscillator module allows to output any number of pulses between 1 and 16,772,215 (16.7 million) to one channel before the next channel becomes active ("bursts"). Adjacent output channels can also be combined to emit pulses at the same time.


The output of each channel can be enabled or disabled. This does, however, not mean that the channel is "eliminated" from the rotary working principle. Instead, the programmed number of pulses is still processed, but simply no signal is present at the output. This is useful to insert time gaps between two bursts.

The synchronization signal


preliminary

Each pulse, that is output at any of the eight channels, can be accompanied by a synchronization signal at the sync out connector of the oscillator module. This synchronization signal can be used to, e.g., indicate proper operation to an amplifier system.

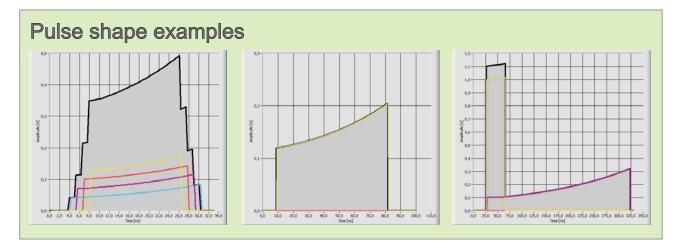

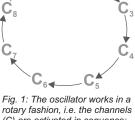

(C) are activated in sequence: 1, 2, 3, 4, 5, 6, 7, 8, 1, 2, 3,...

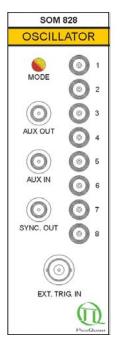

preliminary

preliminary

Specifications

Oscillator module (SOM 828)
Input
Operation modes rotary, programmed sequence of one channel must be completed before next channel is activated, multiple channels can be combined
Oscillator type crystal locked
Base frequencies
Synchronization output
Timing
External trigger input Amplitude
Bursts
Burst length
Waveform generation module (SWM 828)
Input
status control synchronization output Number of waveform generatorsup to 4
Waveform parameters
Time base length
Operation modes
Continuous pulse mode pulse repetition rate from once every minute to 1 MHz and single pulse Burst pulse mode up to 16 million pulses per burst, pulse repetition rate from 196 kHz to 1 MHz, burst repetition rate from once every minute to 500 kHz
Voltage controlled laser module (VCL 828)
Inputs two analog inputs for output from SWM 828, max. superimposed analog voltage: 1 V for full optical power
Optical output
Center wavelength
Output connector FC/APC fiber connector
Output power
Mainframe
Power supply
Computer
Operating system
Please check our website for updated information.
INVISIBLE OR VISIBLE OR VISIB

All Information given here is reliable to our best knowledge. However, no responsibility is assumed for possible inaccuracies or omissions. Specifications and external appearances are subject to change without notice. Trademarks or corporate names are used for explanation and identification, to the owner's benefit and without intent to infringe.

© PicoQuant GmbH, April 2012

Phone +49-(0)30-6392-6929 Telefax +49-(0)30-6392-6561 Email info@picoquant.com

WWW http://www.picoquant.com

preliminary

CLASS 3B LASER PRODUCT IEC / EN 60825-1



PicoQuant GmbH Rudower Chaussee 29 (IGZ) 12489 Berlin Germany

